IKA

designed for scientists

IKA lab dancer

Betriebsanleitung Ursprungssprache	DE	3
Operating instructions	EN	6
Mode d'emploi	FR	9
Instrucciones de maejo	ES	12
Handleiding	NL	15
Istruzioni per l'uso	IT	18
Driftsanvisning	SV	21
Driftsinstruks	DA	24
Driftsvejledning	NO	27
Käyttöohje	FI	30
Instruções de serviço	PT	33
使用说明	ZH	36
取扱説明書	JA	39

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Contents

	Page
EU Declaration of conformity	6
Safety instructions	6
Correct use	7
Unpacking	7
Commissioning	7
Maintenance and cleaning	7
Warranty	7
Technical data	8

EU Declaration of conformity

We declare under our sole responsibility that this product corresponds to the regulations 2014/35/EU, 2014/30/EU and 2011/65/EU and conforms with the standards or standardized documents: EN 61010-1, EN 61010-2-051, EN 61326-1, EN 60529 and EN ISO 12100.

A copy of the complete EU Declaration of conformity can be requested at sales@ika.com.

Safety instructions

In order to be able to use the appliance properly and safely, every user must first read the operating instructions and observe the safety instructions contained therein. Take care of these operating instructions and keep them in a place where they can be accessed by everyone.

The **IKA lab dancer** is a laboratory instrument which conforms to the international safety regulations for laboratory instruments.

Caution: overfilling with liquid may cause the medium to splash out.

You should therefore always use covered test tubes.

The unit is not suitable for manual operation. Observe all pertinent safety instructions and directives, as well as work-safety and accident-prevention regulations for laboratory use.

The appliance must only be used on a non-slip surface.

Grubby support surface diminish the stability of the appliance.

When working with the shaker, the user must select and wear his personal protective equipment according to the mixing hazard category. Defective or inadequate protective equipment can expose the user to the risk of squirting liquids and projectile parts. For this reason, only closed test tubes should be used.

The user may also be at risk if inflammable materials are used in the shaker, or if the trigger contact pressure is too great for glass equipment or the transmitted mechanical shaking power causes the glass to break.

Furthermore, the shaker may only be used to shake materials or material mixtures that the user knows will not react dangerously to the extra energy produced by the shaking. This also applies to extra energy generated by solar radiation during the shaking procedure.

The appliance is maintenance-free and must not be opened.

The **IKA lab dancer** was not designed for use in dangerous atmospheres, for mixing dangerous substances or for use under water.

If the appliance begins to slide, wipe the soft stand feet with a damp cloth.

Correct use

The **IKA lab dancer** is designed for shaking test tubes. Gently push the test tube into the soft attachment to start shaking. Shaking stops as soon as you lift out the test tube.

Unpacking

Please unpack the device carefully and check it for damage. Any damage should be notified immediately to the shipping agent (post office, railway network or logistics company).

As supplied, the package contains:

The **IKA lab dancer**, a plug-in power supply unit and an instruction manual.

Commissioning

Connect the appliance to a socket (Please refer to the information in the "**Technical Data**" section).

Maintenance and cleaning

The **IKA lab dancer** is maintenance-free. It is subject only to the natural wear and tear of components and their statistical failure rate.

Use only cleaning agents which have been approved by **IKA** to clean the device: water containing surfactant / isopropyl alcohol.

- Wear protective gloves while cleaning the devices.
- Electrical devices may not be placed in the cleansing agent for the purpose of cleaning.
- Do not allow moisture to get into the device when cleaning.
- If a different cleaning or decontamination method than the method defined by **IKA** is planned, the user must ascertain with **IKA** that this method does not damage the device.

Warranty

In accordance with **IKA** warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover worn out parts, nor does it apply to faults resulting from improper use, insufficient care or maintenance not carried out in accordance with the instructions in this operating manual.

Technical data

Plug-in power supply unit		
Rated voltage	VAC	100 ... 240
Rated frequency	Hz	50/60
Output voltage	VDC	12
Absorbed power	W	max. 30
Output power	W	7.2
Power input	W	1.5
Drive		
Speed	rpm	2800
Operating time	%	100
Protection class		II
Pollution degree		2
Overvoltage category		II
Protection class acc. to EN 60529		IP 40
Ambient humidity (relative)	%	80
Ambient temperature	°C	+5...+40 (in operation)
Transport temperature	°C	-25 ... +70
Storage temperature	°C	-10 ... +70
Weight	kg	0.55
Permissible shaking weight	kg	0.1
Test tube size (Ø)	mm	30
Dimensions (Ø x H)	mm	100 x 70

Type of movement		orbital
Orbital diameter	mm	4.5
Touch function		yes

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IKA Vietnam Company Limited
IKA India Private Limited